



NMR Tips

VIEWING SPECTRUM DURING ACQUISITION

After at least one block has been acquired

Type **wft aph0** in the command line

Type **lb=10** will add 10 Hz line broadening to the spectrum, redisplay by typing above command

Block size can be set in the acquisition tab prior to starting the acquisition

RESETTING SPECTROMETER WHEN THERE IS A COMMUNICATION PROBLEM

Right-click on desktop and select **open terminal**

Type in **su acqproc** and press return

Press the reset button on the spectrometer control board and wait until all boards reset
(scrolling status lights)

Type in **su acqproc** and press return